

# SAMPLING INSTRUCTIONS

WATER BACTERIOLOGICAL PURITY & NITRATE ANALYSIS

## PAID

DATE	AMOUNT
<input type="checkbox"/> CC	<input type="checkbox"/> CK #
<input type="checkbox"/> CA	


## SAMPLING KIT

Contains two containers, a clear 125-ml plastic bottle for coliform analysis, and / or a 125-ml poly bottle for the nitrate analysis. This kit is designed for coliform and nitrate analysis only. No other tests will be performed from the containers.

Bacteria samples brought in

- **On Fridays**
- **Day before a Holiday**

**Will incur a \$50 surcharge**

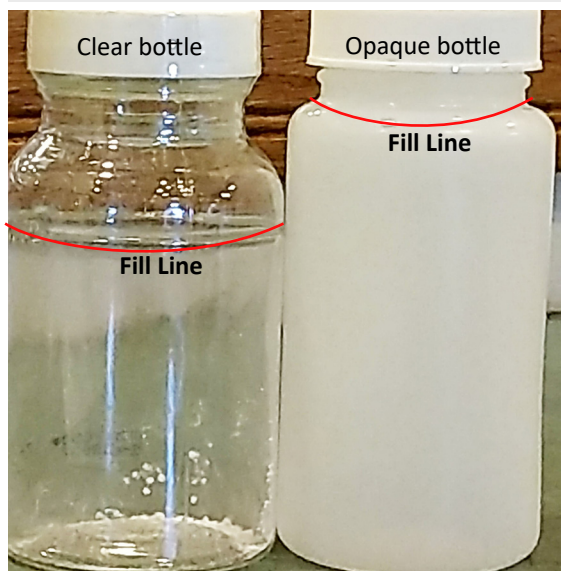

## SAMPLE COLLECTION STEPS

**STEP 1:** Select a cold water tap that supplies untreated or unfiltered water. This ensures a sample that is representative of the source water.

**STEP 2:** Allow the water to run for several minutes until cold.

**STEP 3:** Fill the clear 125ml container to the line and replace the lid.

**STEP 4:** Fill the second opaque container, to the neck and replace the lid.

**Caution: Do not rinse the containers.**

**STEP 5:** Complete the form below and deliver filled containers to Enviro Service, Inc. within **24 hours** of the sampling time.

## FILL OUT

Where sample was collected: \_\_\_\_\_

Date collected: \_\_\_\_\_ Time collected: \_\_\_\_\_

Sampler Name: \_\_\_\_\_

Name: \_\_\_\_\_

Street or P.O. Box: \_\_\_\_\_

City, State Zip: \_\_\_\_\_

Phone: Home \_\_\_\_\_ Cell: \_\_\_\_\_

Work: \_\_\_\_\_ Fax: \_\_\_\_\_

E-mail Address: \_\_\_\_\_

### \*EMAILS:

Will only be sent to you related to business (Including annual test reminders, results, or billing information). You may opt out of e-mails at anytime.

WHITE COPY - LABORATORY    YELLOW COPY - ACCOUNTING    PINK COPY - FILE COPY